

BIOEASY GOAT MILK ADULTERATION RAPID TEST

Bioeasy Goat Milk Adulteration rapid test is a rapid detection for the presence of less than 0.5% of extraneous cow's milk in goat milk sample, not only for fresh goat milk but also for goat milk powder.



RAPID

Results in only 8 minutes



SENSITIVE

Detection of less than 0.5% of extraneous cow milk in any goat milk sample



EASY-TO-USE

No extra devices required, easily perform on site or in lab



SIMPLE

Visually interpretation of results




ACCURATE AND RELIABLE



COST EFFECTIVE

HOW TO USE

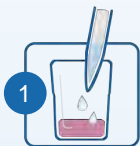



1 Take 200ul sample and mixed



2 Incubate 3 min at 40°



3 Place Test Strip

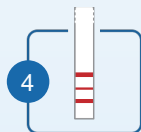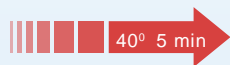

4 Visually Interpretation

RESULTS

Visual interpretation of the result by comparing the intensity of the "TEST" line with the one of the "CONTROL" line.



C—Control Line

T—Test Line

SENSITIVITY

COMPOUND	BIOEASY LOD
Cow milk	0.4-0.6% or 0.1%-0.2%

ORDERING INFO

ORDER CODES	KIT COMPONEN
YREC1001-2	1, Operation instruction 2, 96 test/kit: 12 tubes, each tube with 1 strip of 8 reagent micro-wells and 8 test strips.

STORAGE

At 2-8°C, out of sun light and moisture; Maintained sealed, do not freeze.
Shelf Life: 12 months